



CHROMagar E.Coli 0157

Cat No. LD520

A new chromogenic medium for isolation of *E. coli* 0157 specimens enterohaemorrhagic, allowing direct differentiation of the microorganism.

The medium includes specific chromogenic substrates which allow specific microorganisms become recognizable at a glance by their colony color.

Hy-Laboratories Ltd.
Park Tamar
Rehovot 76326 Israel
Tel : 972-8-9366475
Fax : 972-8-9366474
hylabs@hylabs.co.il
www.hylabs.co.il

Apply procedures identical to those used in classical methods:

1. Direct inoculation from stools on CHROMagar 0157 or subculture after a resuscitation method such as immunocapture on specific beads coated with antibody against 0157.
2. Overnight incubation (24 hours) at 37°C.
3. Rapid and reliable detection and differentiation.
4. Confirm by serological test.

E. coli 0157 - Purple colonies

Most other microorganisms - Inhibited or blue colonies